Synthra Various Radiosynthesizer Product Description and Technical Specifications

Synthra MChelateplus (Catalog No. 013p)

Synthra MChelateplus is a flexible and completely automated radio synthesis system for routine production of a wide variety of Metal- ([⁶⁴Cu], [⁸⁹Zr], [⁴⁴Sc], [⁶⁸Ga], [⁸⁶Y], [¹⁷⁷Lu]) labeled compounds e.g. M-PSMA-, M-DOTA-, M-NOTA-, M-TE2A-, M-TETA-, M-ATSM-, M-PTSM-based radiotracers. Automating the synthesis is simple, with the easy-to-use configuration software SynthraView. The Synthra MChelateplus module offers both, fully automatic and manual modes of operation.

Labeling Possibility Examples

Peptide application: DOTATOC or **DOTATATE** can be labeled by heating 50 nmol of peptide at pH 3.5 – 4.2 for 5 min at 95 °C. For purification, the reaction mixture is passed over a C-18 cartridge to avoid any potential metal breakthrough.

- [⁶⁸Ga] Labeling: The automated synthesis takes about 20 minutes with a decay-corrected yield of about 50 %.
- [¹⁷⁷Lu] Labeling: The automated synthesis takes about 15 minutes with a decay-corrected yield of > 70 %.

[⁸⁹Zr] solution application: Monoclonal antibodies (mAbs) conjugated with DFO can be labeled fully automated at room temperature in 60 min. The reaction mixture can be purified and desalted using a PD-10 SEC column. The automated synthesis takes about 80 minutes with isolated yield of > 70 %.

General Features

- Heating and cooling capabilities
 - One heating zone with cooling capabilities
 - Temperature range: -50 °C 200 °C
- Detectors and controllers
 - Four shielded radiation detectors
 - Two pressure sensors
- ✓ Built-in preparative radio/UV-HPLC system, fixed wavelength LED detector with 255 nm or 280 nm, HPLC column and solid phase extraction (SPE)
- ✓ Self-cleaning system
- Liquid nitrogen trap for radioactive volatiles and to protect built-in vacuum pump



Dispensers and valves


- HR-dispenser (up to 50.000 steps, 2.5/5 mL)
- HPLC pneumatic injection valve (3 mL sample loop)
- Three spare valves for customization
- Chemically inert valves with small dead volume < 35 μL, 5 bar rated
- ✓ Dimensions (w x d x h): 52 × 50 × 48 cm
- ✓ Weight: approx. 35 kg

Synthesis Features

- Eight reagent vials
 - Six small (1 3 mL) and two large (10 15 mL) volume glass vials for reagents
- Three cartridge holders for in-process purification
- SPE unit for final product formulation

Additional Synthesis options

- Product solvent evaporator (Catalog No. 000pse)
- → [¹⁸F] option with [¹⁸O]water recovery system (Catalog No. 013f)

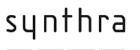

Synthra Various Radiosynthesizer Product Description and Technical Specifications




→ Variable wavelength UV detector (Catalog No. 000vuv)

GMP Features

- ✓ Synthesis files for various tracers
- ✓ GMP compliant. Electronic control and data collection (27/18 channels)
- ✓ 21CFRpart11 & LIMS compatible

Terminal Control

- ✓ A laptop (Win 10 Pro) with preinstalled controlling software SynthraView is included
- ✓ Four digital inputs for communication with external devices upon request

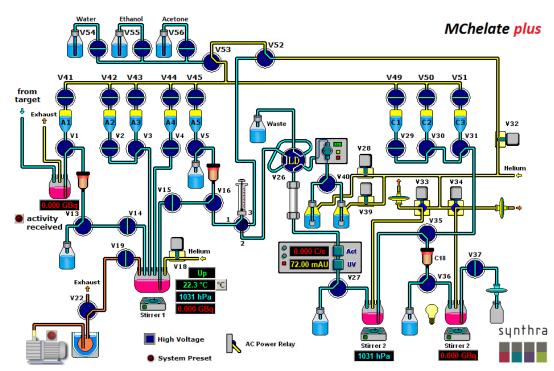

The Graphical User Interface (GUI) of the SynthraView software.

Synthra GmbH reserves the right to modify the information contained herein without prior notice. Copyright © 2022 Printed in the EU

2